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
ANEXO:

CCS175 SERIES OPERATION MANUAL
THORLABS SPECTROMETER CCS175

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THORLABS
Spectrometer
CCS175
 $\lambda = 500 - 1000 \text{ nm}$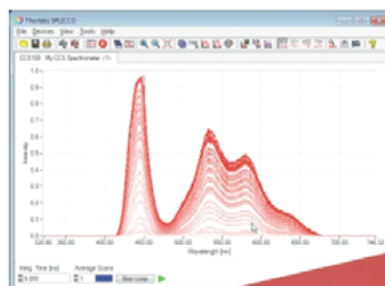

1. GENERAL INFORMATION

The CCS Spectrometer Series is designed for general laboratory use. Integrated routines allows averaging, smoothing, peak indexing, as well as saving and recalling data sets. The initial setup is simple to complete. Following installation of the software, the CSS spectrometer is ready to use. Simply plug it into a USB 2.0 port and run the application software SPLICCO. The remainder of this manual is devoted to the setup procedure and features of the fiber spectrometer. A troubleshooting section and detailed specifications of the various components are provided to further assist. The description of the instrument driver commands can be found in the VXIpnv VISA instrument driver package.

Application software SPLICCO

SPLICCO is an acronym for “**SP**ectrometer and **L**ine **C**amera **C**ontrol”. This software can be used for acquiring direct, transmittance and absorbance measurements in conjunction with Thorlabs line cameras and spectrometers.

After the installation the software is able to communicate with all Thorlabs CCS spectrometers. Additionally, two virtual devices are included: a line camera and a spectrometer, to demonstrate the functionality of SPLICCO.

2. CCS SERIES OPERATING INSTRUCTIONS

Start the GUI



To start SPLICCO click on the desktop icon  or select 'Programs' via the START button in the Windows task bar and navigate to 'All Programs / Thorlabs / SPLICCO / SPLICCO'.

CCS Spectrometer with all user relevant ports and signal LEDs


Connecting a device

1. To start a measurement with a CCS spectrometer, connect it to a USB port of your PC with the supplied cable.
2. Select 'Connect...' from the Devices menu or click to the Connect icon  from the tool bar.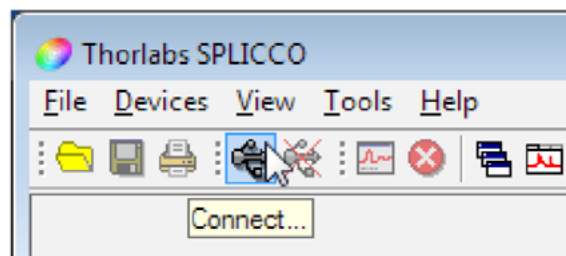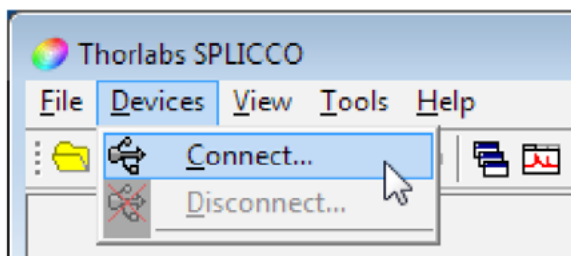

The following window appears and shows all connected devices and additionally two virtual devices. Now you can select a device to be used. A panel will be created according to your selection by default. If the “Create a device window automatically” option is not checked, please use the according panel icon from the main interface.

Press “Cancel” to leave this dialog and “Rescan” to scan the system again for new devices.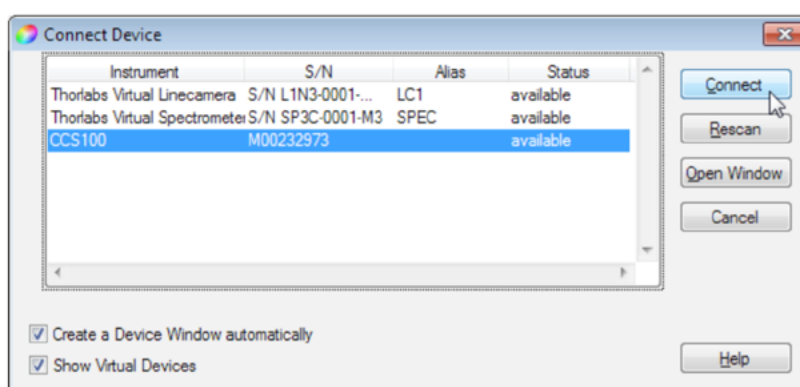

Every device can only be opened once. Devices already opened by SPLICCO are marked with the “running” status and are grayed out in the device selection dialog. Devices used by another application than SPLICCO are marked with the “locked” status.

Press “Open Window” to switch to the “Open Window” dialog to connect a window to an already running device.

Furthermore, you can start a virtual spectrometer, which can simulate a spectrum. Through this feature you can familiarize yourself with SPLICCO, without the need of a light source or signal. You can select to display or hide those virtual devices by checking or unchecking the “Show virtual devices” box.

For a detailed description about the virtual devices refer to section Virtual devices.

Integration time

The integration time represents how long the CCD interacts with incoming light. CCD pixels act like light buckets, gathering photons. The integration time displays the duration for which the bucket is open. For very bright sources, low integration times are required, whereas for weak sources, longer integration times should be used. As in the light bucket analogy, CCD pixels can be overfilled (“blooming”). This is called saturation and will cause the output to be misleading.


If no intensities are displayed, please enter the shortest integration time and increase it continuously until an intensity curve is displayed. As mentioned above, CCDs are very sensitive, and if over-exposure occurs, no intensity can be displayed.

Also, please make sure the background correction is disabled (see section Background Correction)

Integration time can be set via the control on the lower left corner of the device window. The supported range is defined by the CCS Spectrometer and ranges from 10 μ s to 60 seconds. The integration time input window uses milliseconds, therefore the values of 0.01 ms – 60000 ms have to be used to cover the range. A window, which only shows a loaded from a file spectrum does not offer those controls in the left bottom corner.


A change of the integration time affects all windows connected to this device, which are then updated to show the same integration time.

Higher integration times results in higher peaks in the measurement data.

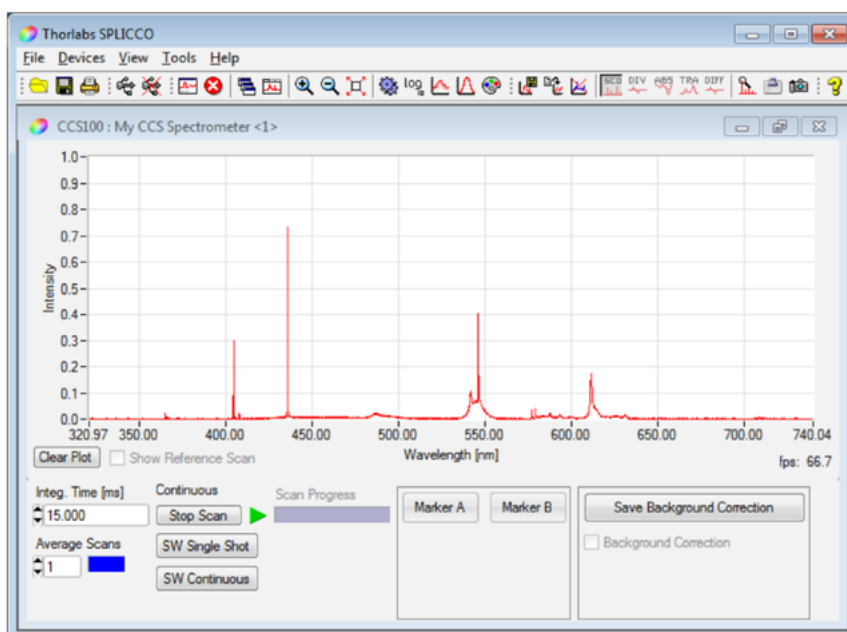Loading data


Previously captured and saved measurement scans can be reopened with SPLICCO without connecting a device. This can be done from the File menu by selecting “Open...” (📁 icon) or “Import CSV...”.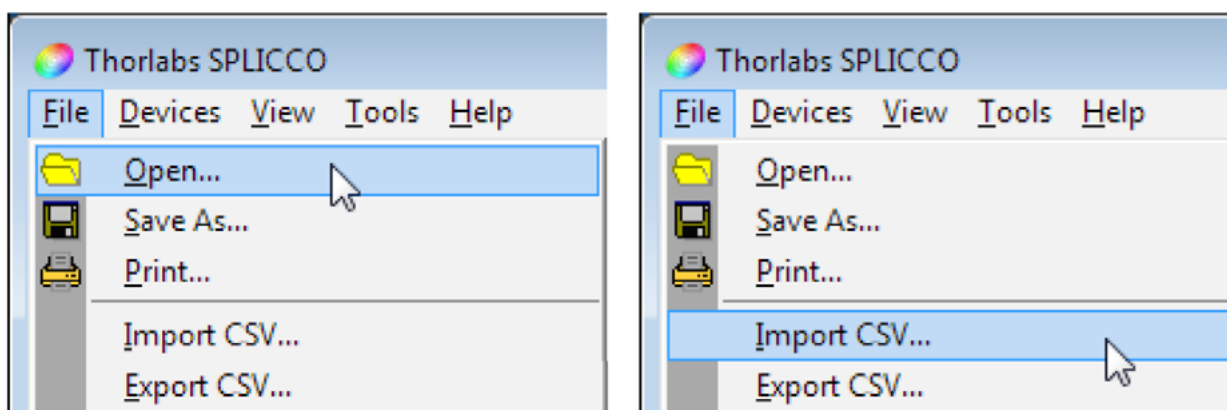

Opening a file

SPLICCO can load most JCAMP-DX files with file extension *.jdx either as a reference to use with the live measurement data or in an individual window to show formerly saved data.

To open a *.jdx file, choose “Open” or click to the  icon. Choose the appropriate file using the Browse button, the selected file is shown in the preview window. Eventually, saved comments are shown in the “Comments” field.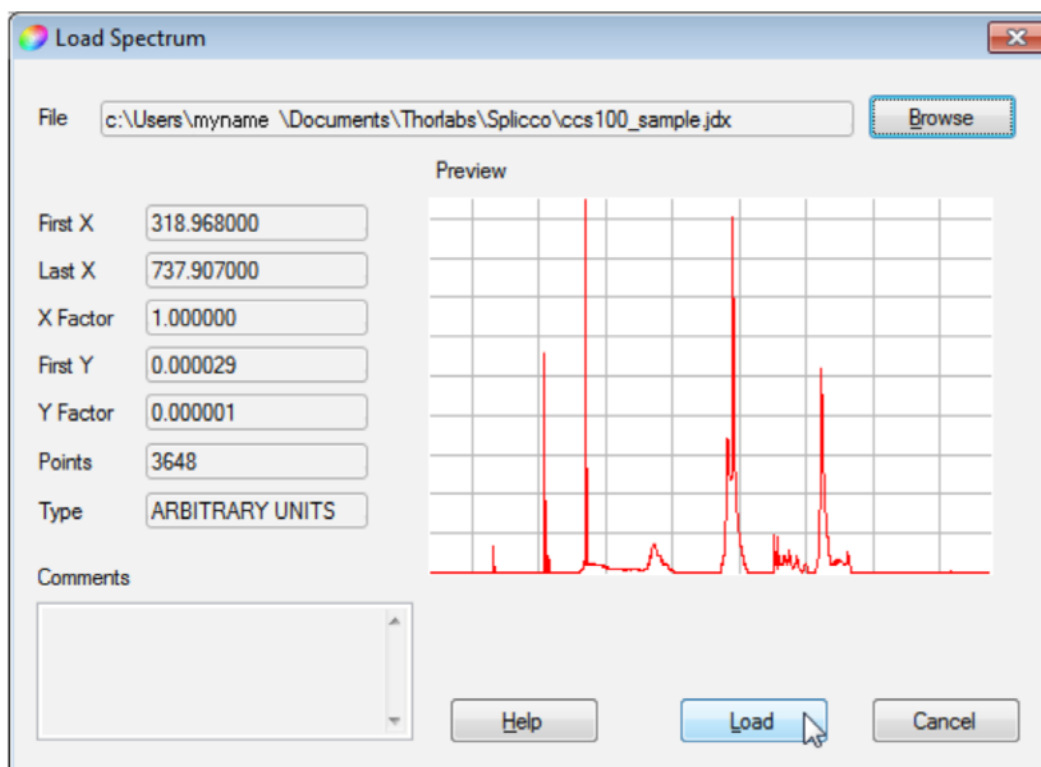


X and **Y** Factors represent the resolution. The resolution of the X axis is 1 pixel, the resolution of Y axis results from the resolution when the file was saved. By pressing Load the curve is opened in a new window.


The **Load** function is used also for loading a reference scan to the current live window. Details are explained in section References.

Import data


To import measurement data from a *.csv file select ‘File -> Import CSV...’ from the menu.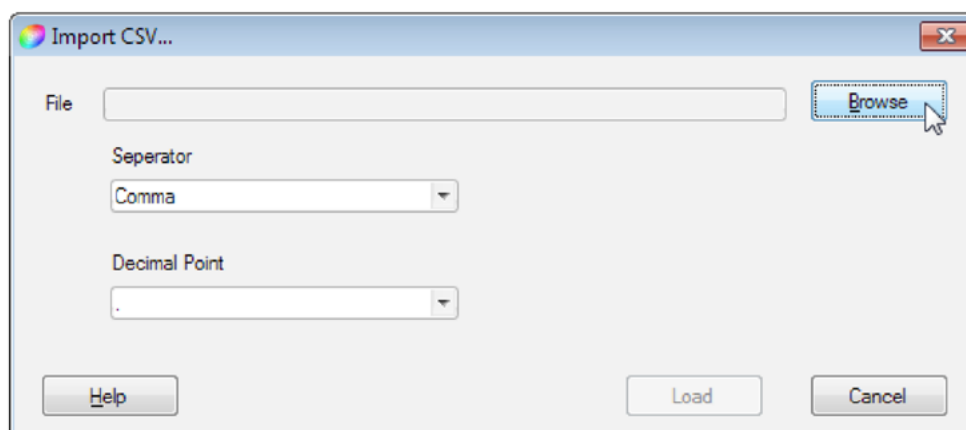

Please specify the character separating the x & y columns and choose which character marks the decimal position in the appearing window. Then click to **Browse** to select the required file:

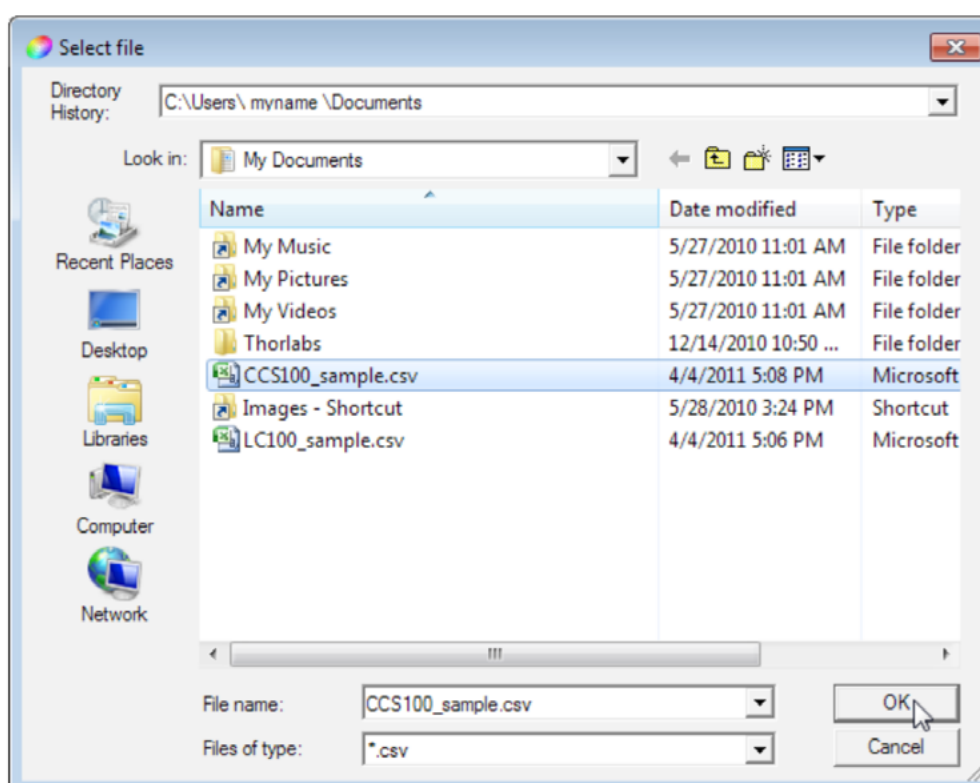Click OK to confirm:


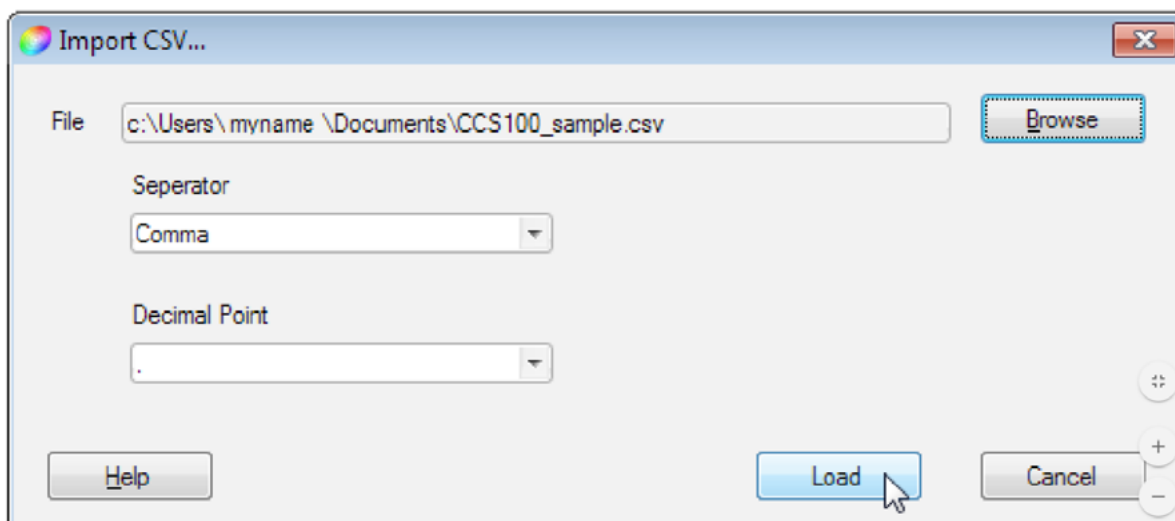
Clicking to **Load** imports the data to a new window.

Saving data

SPLICCO can save data either in JCAMP-DX or CSV format.

To save the measurement data to a JCAMP-DX file select 'Save As ...' from the **File** menu or click the button  from the toolbar. A file dialog window appears and you can choose the filename and directory.

The file extension of this file is *.jdx. Additionally, you can choose the range to store and the resolution, which defines how many values are interpolated and saved to file. "High" means factor 2, "Low" means factor 0.5 and "Average" means factor 1. Text entered in the "Comments" field will be stored together with the data.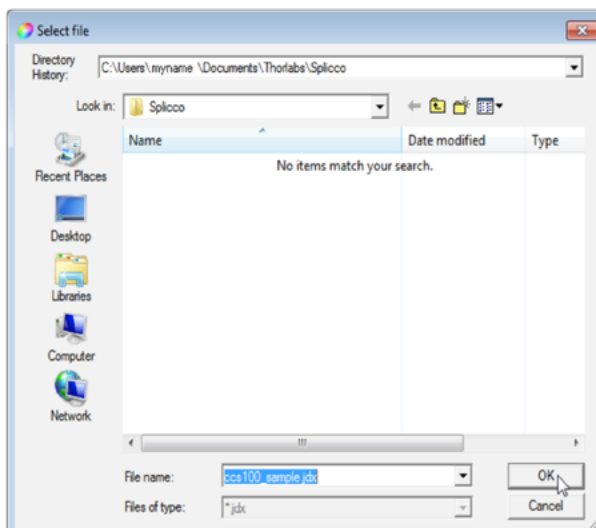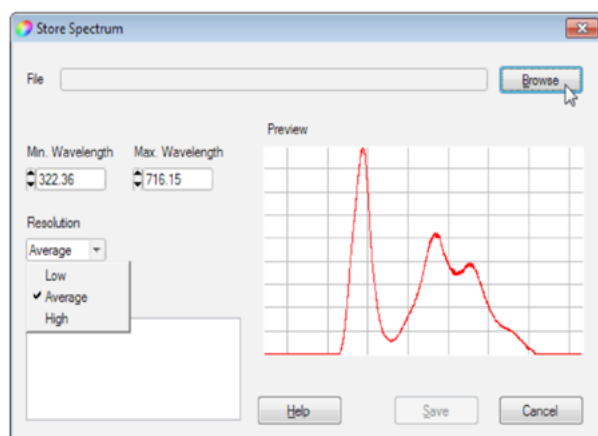


Export Data

Measurement data can be exported to a *.csv file for use with e.g. Microsoft Excel™ or MathLab™. To export the current measurement data to a *.csv file select ‘Export CSV...’ from the **File** menu. A popup panel appears to choose the target directory, filename and the characters for “**Separator**” and “**Decimal Point**”. Reference data can be handled in the same way.

Device Settings

The Device settings dialog can be opened from the **Device** menu.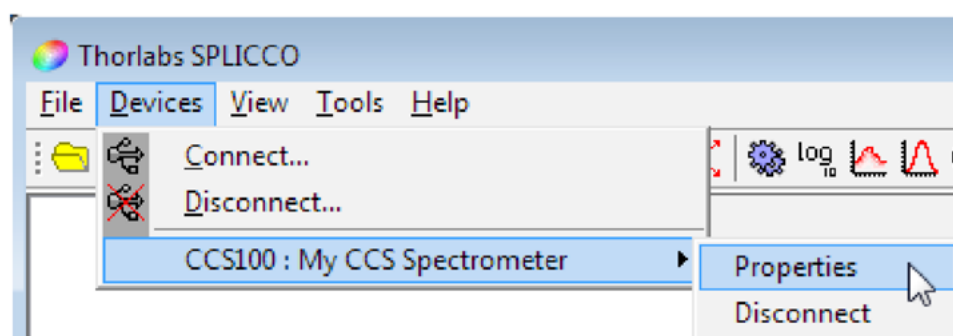

Device Info

The tab Info contains information about manufacturer, device name, serial number and firmware revision.

Trigger mode

SPLICCO is able to generate internal trigger signals (SW) or use external trigger signals to take readings at defined time intervals.


You can set the trigger mode in the **Device Settings**, tab 'Common'.

The control offers five trigger modes: SW Trigger Continuous, SW Trigger Single Shot, External Trigger Continuous, External Trigger Single Shot and the Idle mode. According to the trigger mode the status symbol and the trigger button in the bottom of each device windows changes. The trigger buttons labels shows the possible option, e.g. "Stop Loop", "Scan 1x", "Arm Trigger" or "---".

Averaging

Very noisy or weak signals can be amplified by adding several scans, which is known as averaging. SPLICCO provides two kinds of averaging - Gliding Average and Block Average. The averaging mode can be set in the Device Settings panel in the 'Common' tab.

The number of scans to average can be set in the bottom of the active panel:

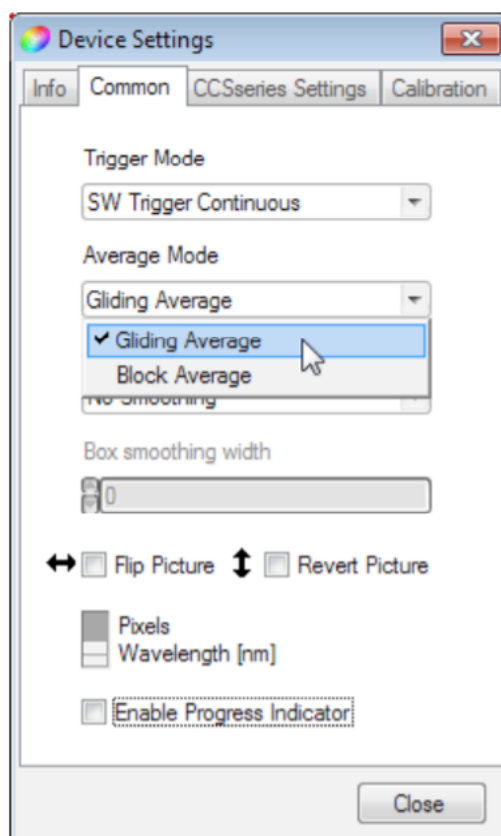To the right of the number of scans to be averaged a status box is displayed, indicating the fill level of the buffer used for averaging.

Gliding Average

This method averages over the most recent number of scans and is being updated with every new scan. The advantage is that the graph is being updated with every scan.

Example: The number of scans to be averaged is set to 10. After starting acquisition, the software calculates the average out of the first two data sets, then out of the first three sets and so on until the desired number (10) is reached. Then the first data set will be subtracted and the newest data set will be added to calculation of average. This can be seen also in the buffer fill level - it grows up to the max and stays there.


Block Average

This method accumulates a number of scans, after that calculates the average, displays it and starts the averaging process from beginning. The display is updated only after n scans.

Display mode

CCS100 Series Spectrometers can use an internal wavelength calibration file to show the range in nanometers instead of pixels. SPLICCO supports two display modes regarding the scaling of the x axis - the x axis is shown in either nano-meter or pixel.

Click to  icon in the toolbar and open the 'Common' tab: flip the switch to change between the display modes.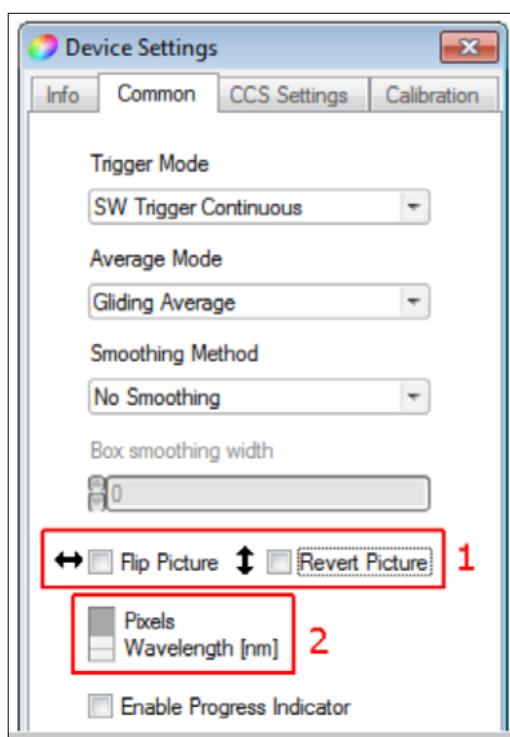


Cursor control

SPLICCO provides a cursor for instantaneous readout of wavelength and amplitude (intensity). In order to use this feature, right click on a graph and choose “Cursor mode”. The “Cursor mode” disables the “Zoom mode” and a vertical line appears. With the first left click on the graph a small window appears right beside the mouse cursor and shows the actual x-value and y-value. The precision is always the same as the precision of the corresponding axis.


Peak finder

SPLICCO offers the possibility to find peaks in an actual measurement. The peak finder can be selected by right clicking on the actual window and selecting “Peak finder...”


In order to specify the relevant peaks, a peak width and a peak threshold can be set. A higher threshold and a higher peak width will reduce the found peaks. By pressing the “Calculate” button the peaks specified by those two parameters are calculated and the list is filled with the found peaks.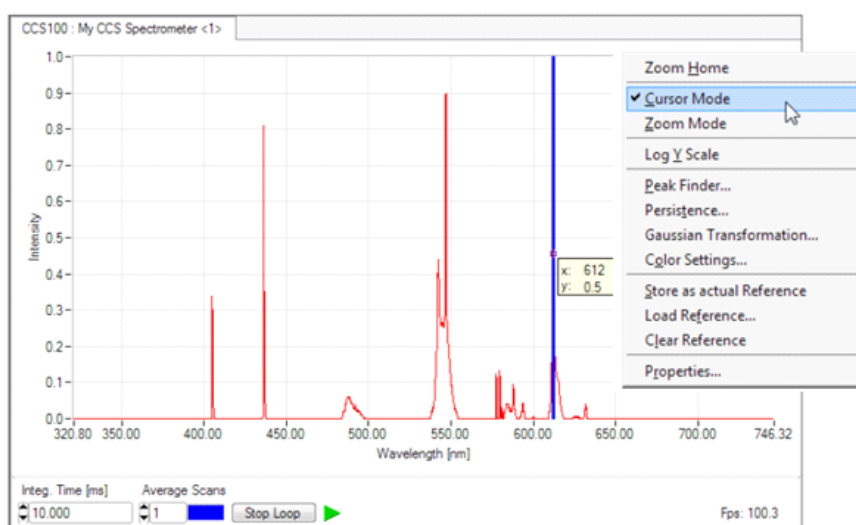

Cursor control

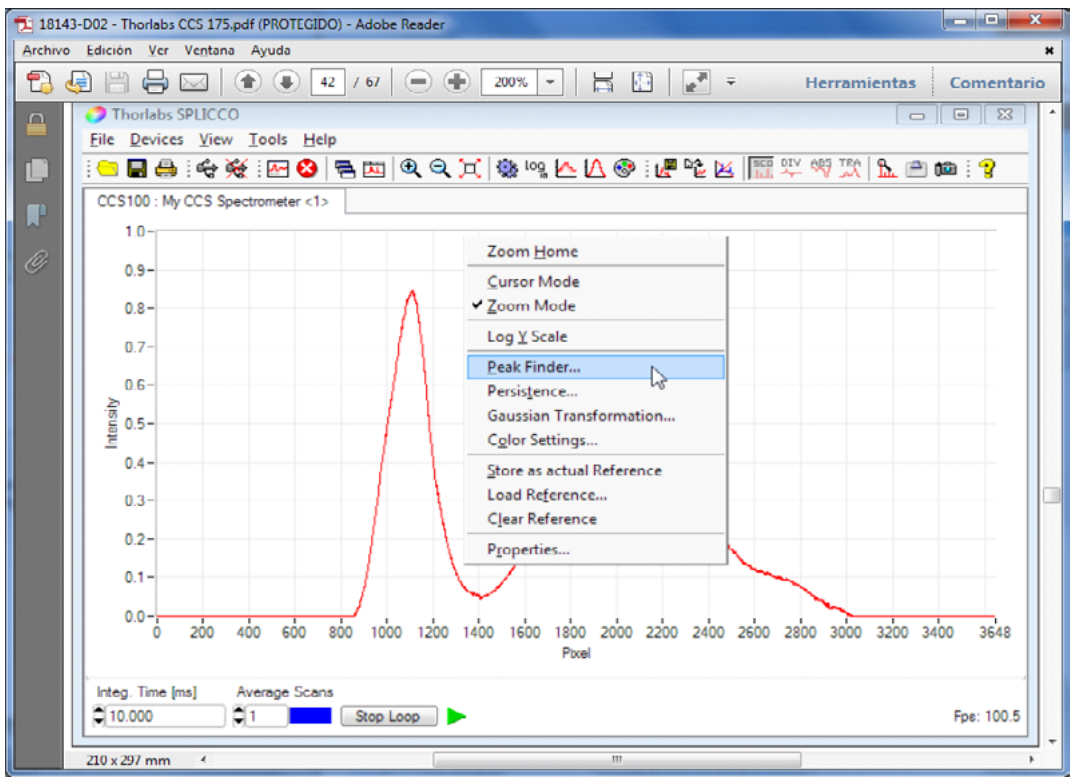Peak finder

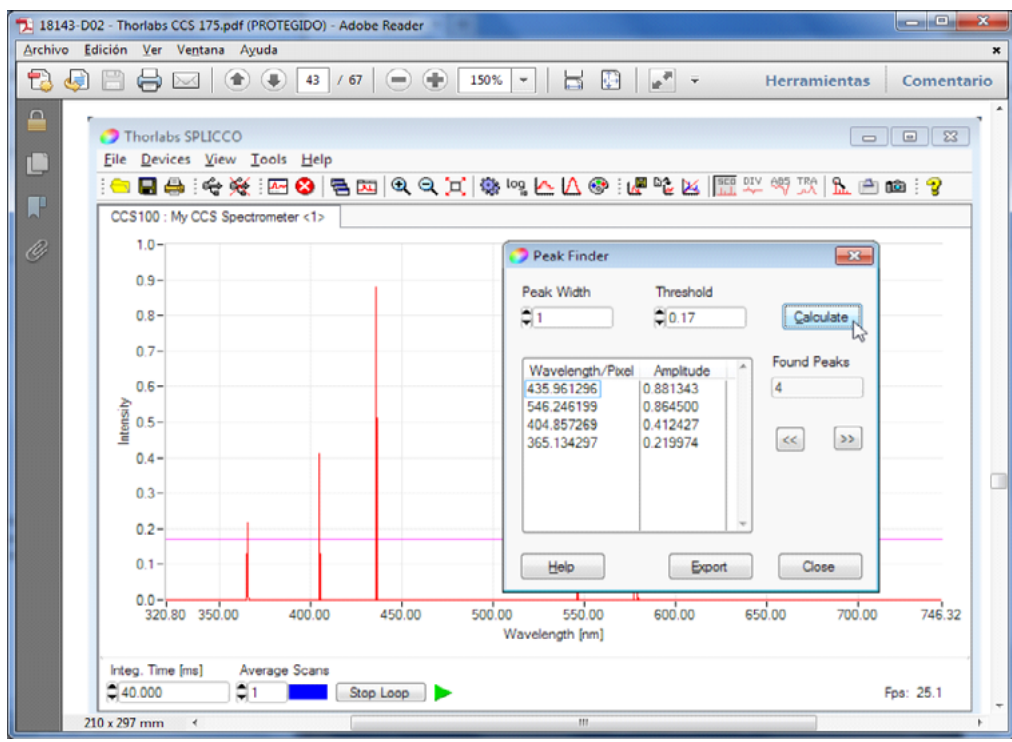Specify the relevant peaks

The found peaks can be sorted by their location (“Wavelength/Pixel”) or by intensity (“Amplitude”) - just click to the desired header - and can be iterated with the help of the arrow buttons (“<<” and “>>”).

Only one peak can be marked at the same time, but you can export the full list into a tab separated text file with the “Export” button.

Logarithmic Y Scale

The intensity axis can be scaled linearly or logarithmical. Default display is linear, in order to switch to logarithmic just click the icon in the toolbar. Alternatively, the **Log Y Scale** can be activated via the **View** menu or from the dialog after right clicking to the diagram area.

References


An actual scan can be stored as reference. This reference will appear in a different color (see Color Setup) and remains unchanged during the current SPLICCO session unless cleared.

Click the  icon in the toolbar, or select ‘**Store as actual Reference**’ from either the **Tools** menu or from dialog, which appears after a right click on the diagram area.


Delete a reference by selecting ‘**Clear Reference**’ or by clicking the  icon. A scan can be saved as a *.jdx file for later use.

Save a reference

- » Store the actual scan as reference. (: In order to distinguish the reference scan in the following screenshot, the integration time was changed after defining the reference. Actually, the actual scan and the reference are congruent).

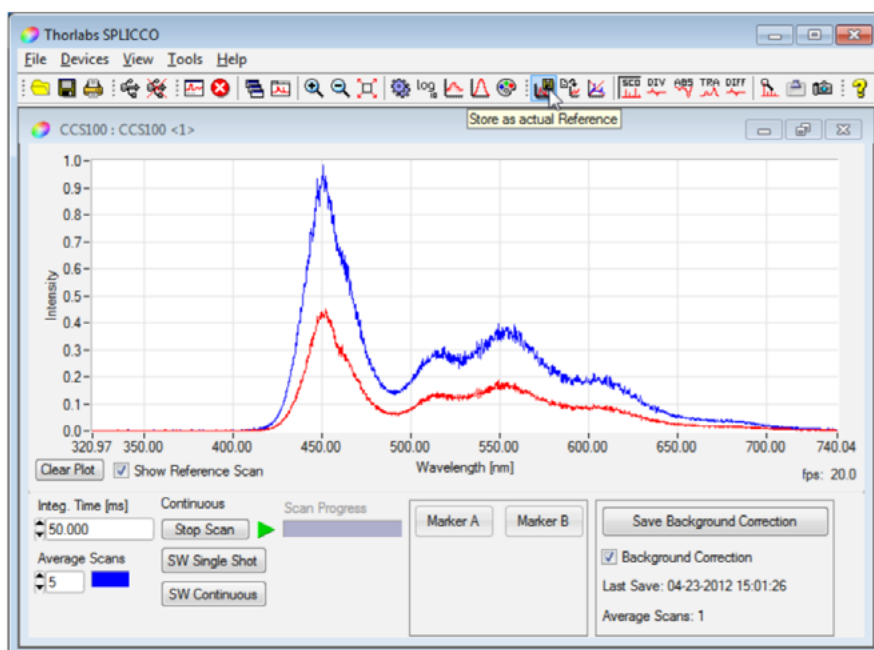
- » Open the 'Save As...' menu, choose a location and file name, select '**Data to store: Reference**' and click **Save**.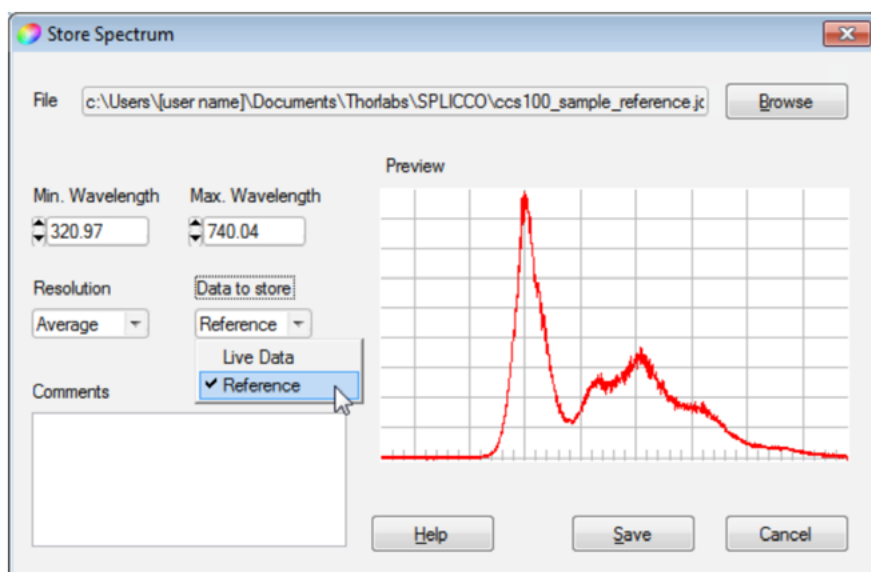

Load a reference

- » Click the  icon in the toolbar, or select 'Load Reference...' from either the **Tools** menu or from dialog, which appears after right clicking the diagram area. A dialog box opens:

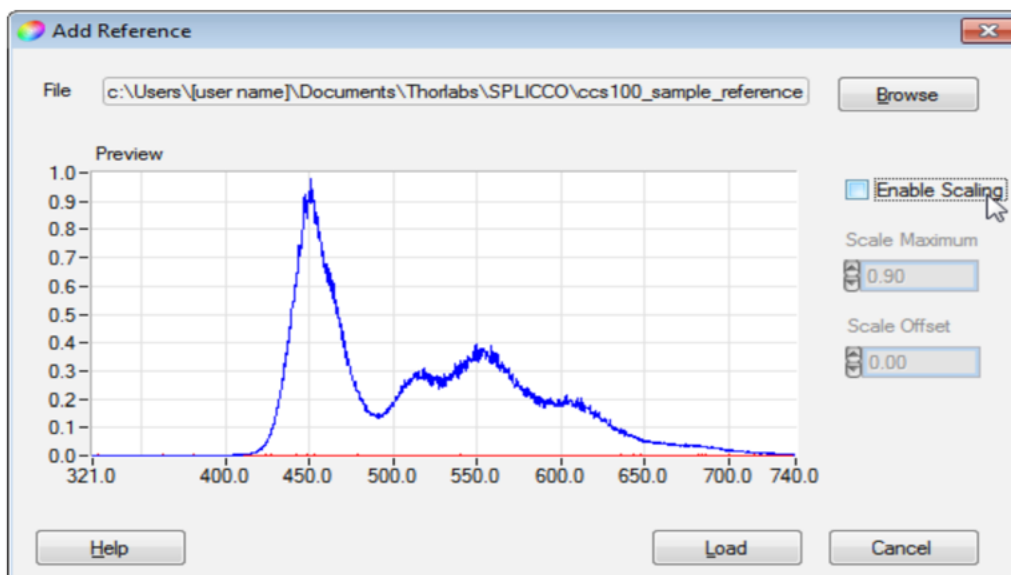- » Browse for the file location, select the desired reference file and click **Load**.
- » The reference curve is copied into the actual window and can be used for future calculations.

Calculations with references

SPLICCO offers the possibility to recalculate the live measurement with a reference curve. This reference curve can be a sampled live curve or a curve loaded from a JCAMP-DX file as long as the ranges are compatible to each other. It is possible to calculate with a reference curve that only partly matches the range of the live measurement. In this case only the matching range is calculated and shown.

There are four modes available:

Scope: The normal view

Relative difference: The relative changes with respect to reference or background

Absorbance: The absorbance view

Transmittance: The transmittance view

An example shall illustrate that. As reference, a spectrum in the range from 420 to 660nm of a Thorlabs MCWHL2 mounted cold white LED was taken. Then, into the light path a Thorlabs NE06A neutral density filter (optical density 0.6, transmission 25%) was inserted.

Scope view

The scope view is the standard view in SPLICCO. The reference curve is shown as well as the original measurement data, no calculations made.

The blue curve is the reference (LED spectrum), the red curve represents the spectrum of the LED with the ND filter.

Relative difference measurement:

The relative difference view shows for each wavelength the live measurement values (I_{meas}) divided by the reference curve values (I_{ref}), so that only changes in the spectrum are displayed. The original reference curve is still shown to allow an easy comparison. The original reference curve is still shown to allow an easy comparison.

Transmittance:


The transmittance view shows the light transmission by a sample in %: Please make sure, the Y axis is scaled to a convenient for display value (here to 100%). The original reference curve is still shown 28 for comparison.

Absorbance:


The absorbance A , also known as optical density OD , describes the light absorption by a sample:


The original reference curve is still shown to allow an easy overview.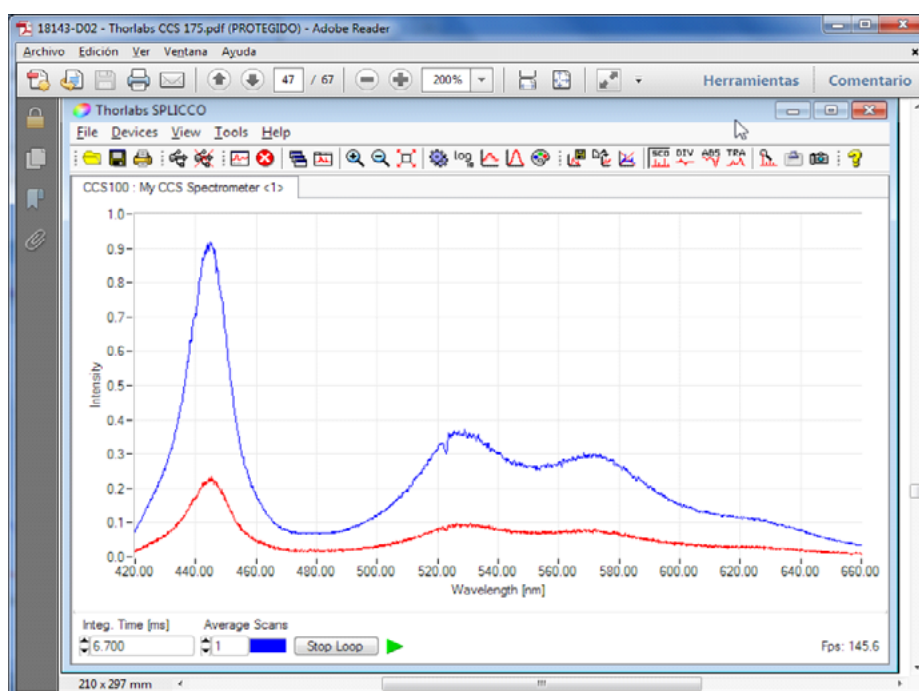


Scope view

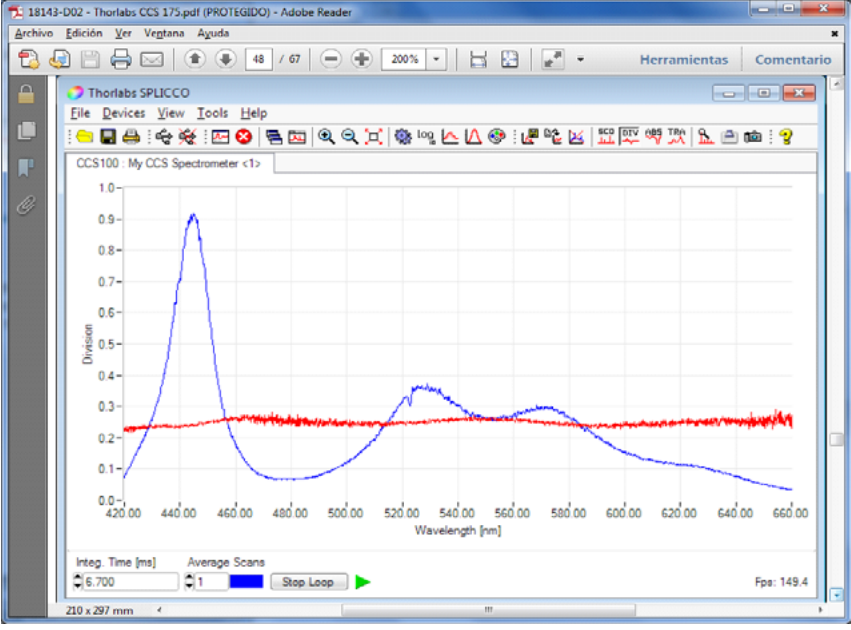Relative difference measurement

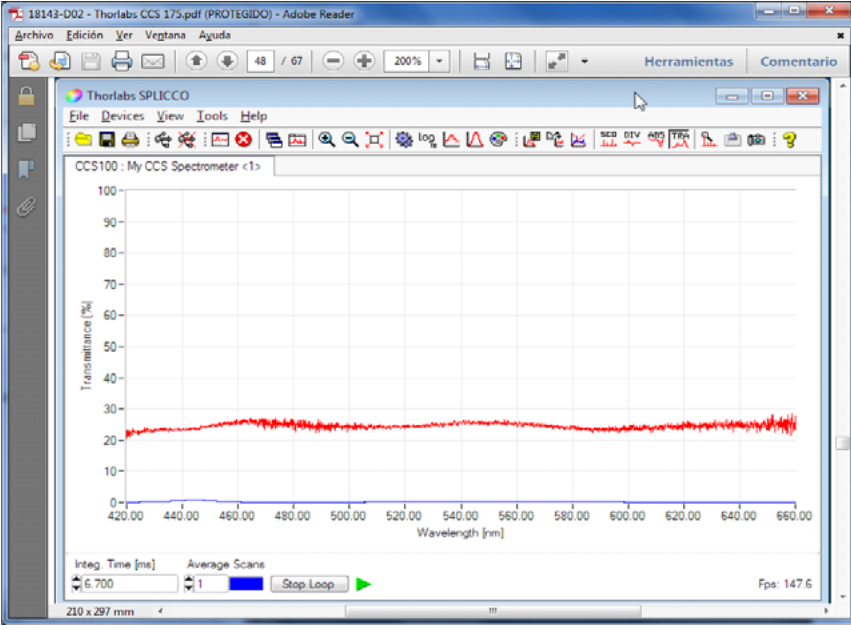Transmittance

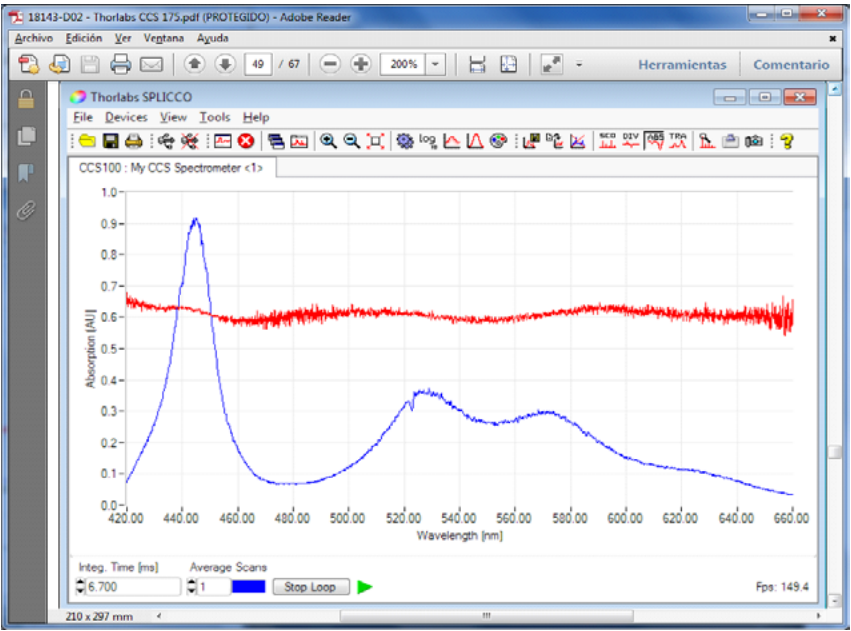Absorbance

Copy to Clipboard

The actual scan data can be copied to the clipboard via the menu **Tools**, choose **Copy to Clipboard**.

These data can be pasted to a different application, e.g. Microsoft EXCEL[®], for further processing. The data are comma separated pairs of values, where the first value represents the pixel number (wavelength in nm), the second - the intensity measured at this pixel.

Snapshot

A **Snapshot** of the actual scan includes the actual scan diagram area and a header:

- » device type and device label
- » date and time
- » user name
- » settings for integration time and averaging counts

To make a snapshot, choose **Take a Snapshot** from the menu **Tools**. A Select File dialog appears, asking for a file name and format. Snapshots can be saved as *.bmp, *.jpg or *.png files.


3. APPLICATION NOTE

This section hopes to aid the user, by giving tips or hints on how to obtain the best measurements with the CCS spectrometers. It begins with a few general suggestions and finishes by giving a choice of question with solutions.

To obtain a good spectrum with minimum noise the light reaching the spectrometer should be maximized. Therefore, a clean fiber with minimum bends is of high importance. Fiber collimation packages, which physically fit to the SMA fiber, but are not designed for Multimode fibers might be less efficient than just the SMA connector in front of the sample.


In some cases a small variation in the diameter of the SMA ferrule, might lead to a tilted fixture of the connector in the fiber port. In this case loosen the fiber connector a few turns and reconnect it.

For a good signal to noise (S/N) ratio, it is recommended to have the spectral data of interest in between 70% and 95% of the intensity scale (0.7 - 0.95). This ensures that the signal is above the background noise, but not overexposed. You can identify an overexposed signal by the increased line width and a characteristic plateau of the signal maximum.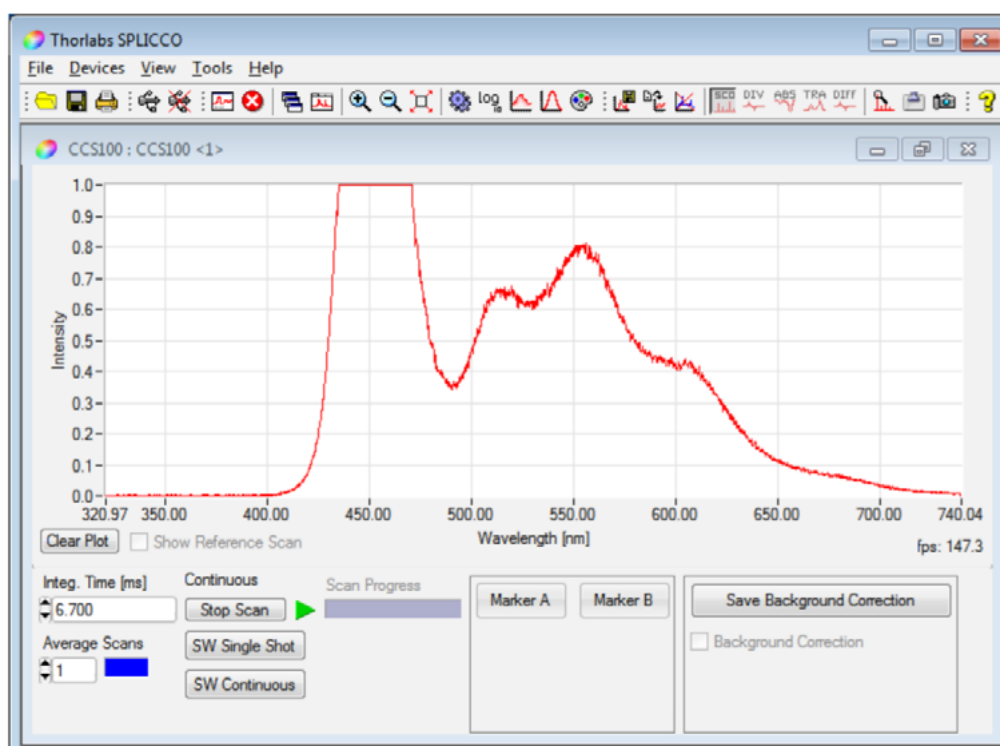


If your signal looks similar to the sample in the above figure, decrease the integration time to get a maximum peak intensity in between 0.9 and 1.0. If your signal is still overexposed at the minimum integration time of $10\ \mu\text{s}$, it is recommended to use neutral density filters or other optical attenuator in front of the light source to decrease its intensity.

If the signal you are evaluating is very small, you can increase the integration time. As shown in the next figures, this way the peak intensity raises from ~ 0.018 (equal to 1.8% of the max. intensity) to about 0.7 (70%), which is a recommended value.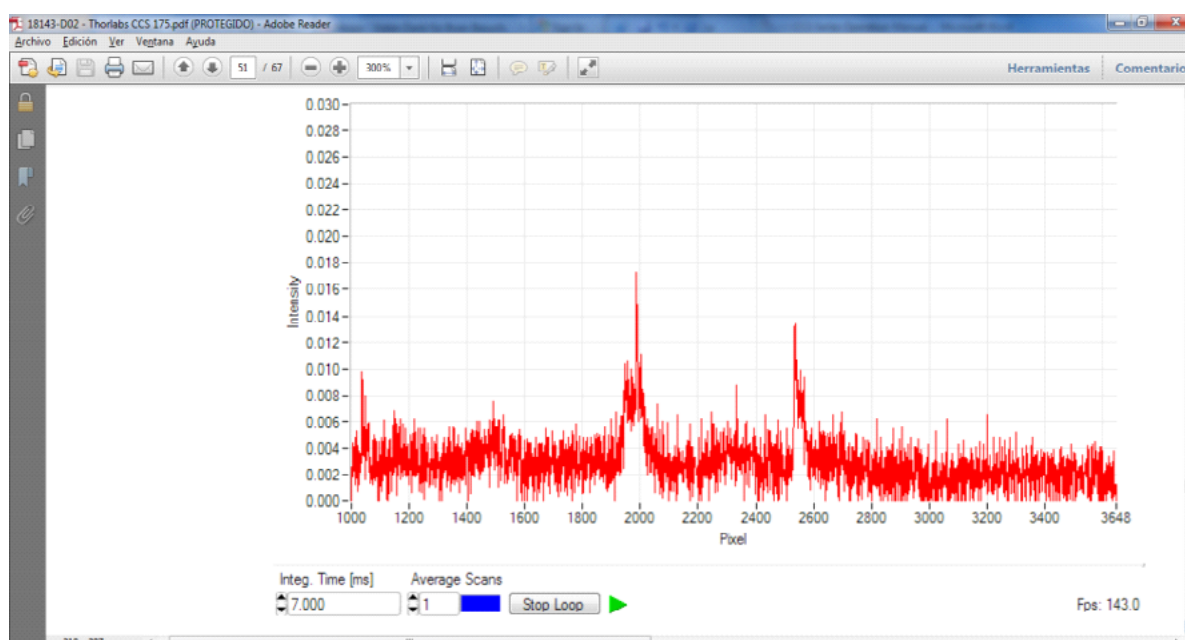

Scan, recorded at 300ms integration time

If you are already at the maximum or if other parameters of your experiment do not permit long exposures, you can also use the averaging function to raise your signal above the noise. The result can be seen in the following figures. Please note, that the intensity is around 0.018 which corresponds to 1.8% of the available intensity scale.

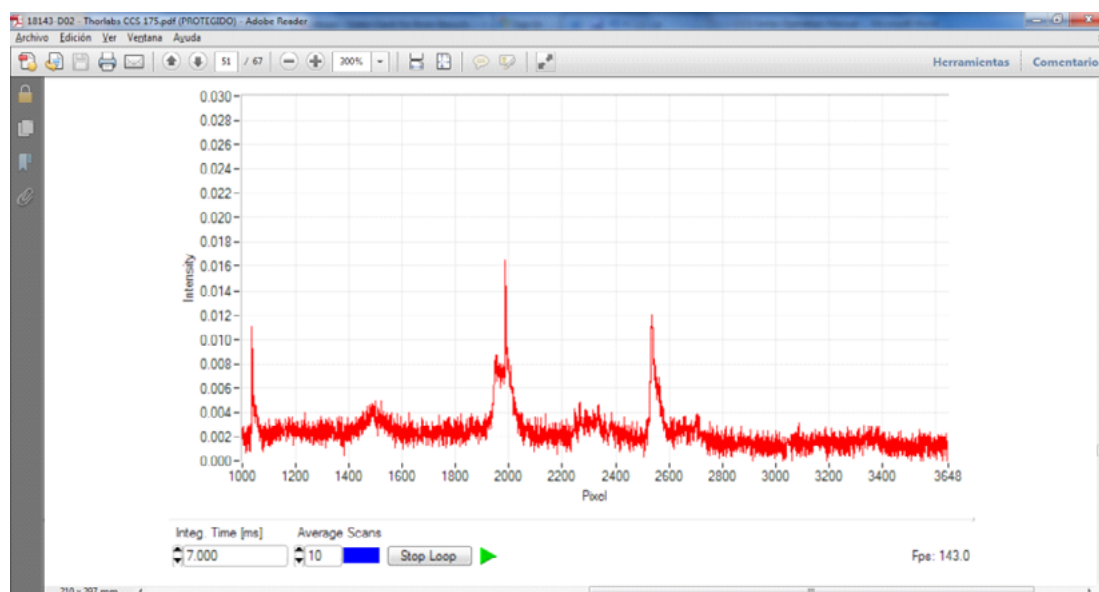Scan, recorded at 7ms integration time and 10 averages

Blooming

A property owned by all CCD sensors, as used in the CCS spectrometers, is blooming.

Strongly overexposed pixels tend to discharge neighboring/adjacent pixels, even if they are not illuminated. This can be seen in the spectra by an increased line width, which could lead to misinterpretations of the signal. You can avoid this effect by decreasing the integration time or using optical attenuators, as described above.

Maintenance and Service

Protect the CCS spectrometer from adverse weather conditions. The CCS spectrometer series is not water resistant.

Attention

To avoid damage to the spectrometer, do not expose it to spray, liquids or solvents!

The unit does not need a regular maintenance by the user.

If necessary the unit and the display can be cleaned with a cloth dampened with water.

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D.R. © 2020, Universidad Nacional Autónoma de México, Avenida